


User Manual



# PH180

## Waterproof pH Meter

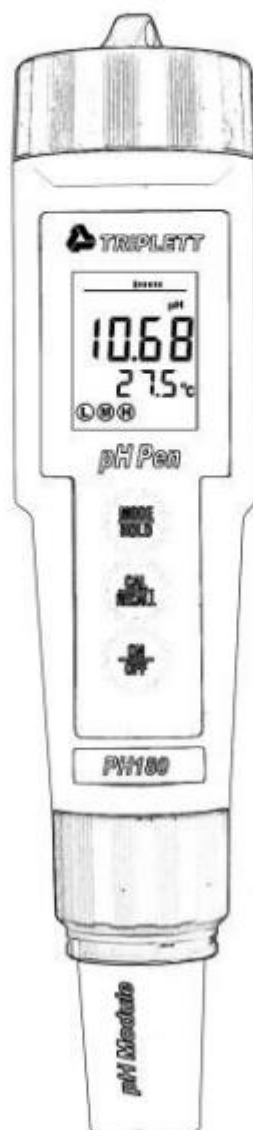

20120067

**DIREKTRONIK**

Kontakta Sverige:  
[www.direktronik.se](http://www.direktronik.se)  
[info@direktronik.se](mailto:info@direktronik.se)  
Ring: 08 52 400 700

Kontakt Norge:  
[www.direktronik.no](http://www.direktronik.no)  
[info@direktronik.no](mailto:info@direktronik.no)  
Ring: 55 50 91 50

# Introduction

---

Thank you for selecting the Triplett PH180 pH meter. This instrument is designed for high accuracy pH testing for solids, semi-solids and liquids. This device is shipped fully tested and calibrated, and with proper use, will provide years of reliable service.

## Features

---

- Measures from 0 to 14 pH with 0.01pH resolution
- ATC (Automatic Temperature Control) compensates for temperature changes
- Dual display for pH and Temperature
- Analog bargraph for viewing trends
- 1-, 2-, or 3- point calibration
- RENEW warning when probe needs cleaning or replacement
- Waterproof (IP57)
- Auto power off (after 10 minutes) with disable
- Includes meter with flat surface pH probe and protective cap, four(4) CR2032 batteries, cup with cap and neckstrap

***DIREKTRONIK***


**Kontakta Sverige:**  
[www.direktronik.se](http://www.direktronik.se)  
[info@direktronik.se](mailto:info@direktronik.se)  
Ring: 08 52 400 700

**Kontakt Norge:**  
[www.direktronik.no](http://www.direktronik.no)  
[info@direktronik.no](mailto:info@direktronik.no)  
Ring: 55 50 91 50

# Description

## METER DESCRIPTION

1. Battery compartment cap
2. LCD Display
3. MODE / HOLD button
4. CAL / RECALL button
5. ON/OFF button
6. Electrode collar
7. Electrode

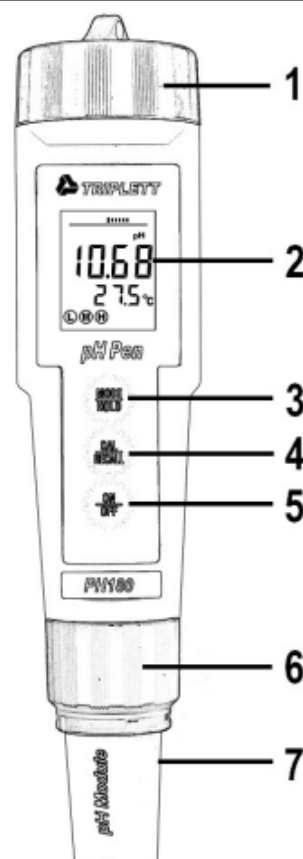

## LCD DISPLAY

1. Bar graph reading
2. Measurement Units
3. Main Display
4. RENEW Indicator
5. HOLD Indicator
6. RANGE Calibration Indicator
7. Low Battery Indicator
8. Temperature Display



## CAUTION

This device is not a toy and must not reach children's hands. It contains hazardous objects as well as small parts that the children could swallow. In case a child swallows any of them, please contact a physician immediately

Do not leave batteries and packing material lying around unattended; they can be dangerous for children if they use them as toys

In case the device is going to be unused for an extended period of time, remove the batteries to prevent them from draining

Expired or damaged batteries can cause cauterization on contact with the skin. Always, use suitable hand gloves in such cases

See that the batteries are not short-circuited. Do not throw batteries into the fire.

## Overview

---

### pH Overview

pH is a unit of measure (ranging from 0 to 14pH) indicating the degree of acidity or alkalinity of a solution. pH tests are the most commonly performed measurements in water analysis and reports the negative log of the hydrogen ion activity of a solution which is an indicator of acidity or alkalinity. Solutions with a pH less than 7 are considered acidic, solutions with a pH higher than 7 are known as bases, and solutions with a pH of exactly 7 are neutral.

The pH scale is logarithmic so, for example, if sample A is 1 pH less than Sample B, this means that Sample A is 10 times more acidic than Sample B. A difference of 1 pH represents a ten-fold difference in acidity.



## Getting Started

For new meters, remove the battery cap and then remove the battery insulating strip.

Remove the cap from the bottom of the PH180 to expose the electrode glass surface and reference junction

Before first use or after extended storage, soak the electrode (with its cap removed) in a pH 4 solution for about 10 minutes

White KCL crystals may be present in the cap. These crystals will dissolve in the soak or they can be simply rinsed with tap water

Always calibrate close to the expected measurement value

A sponge is located in the electrode protective cap. Keep this sponge soaked with a pH 4 solution to preserve Electrode life during storage

## Replacing Electrodes


The PH180 is shipped with an electrode attached. Electrode life is limited and is dependent on (among other factors) frequency of use and care. If the electrode needs to be replaced, follow these steps for removing and connecting electrodes.

1. To remove an electrode, unscrew and completely remove the electrode retaining collar.
2. Gently rock the electrode from side to side, pulling it away from the meter, until it disconnects.
3. To attach an electrode, carefully plug the electrode into the meter socket (note that the electrode connector is keyed, ensuring proper connection).
4. Secure the electrode in place by tightly turning the collar in place. (a rubber gasket seals the electrode with the meter).

## Automatic Electrode Recognition

When the PH180 is turned on, it recognizes the type of electrode that is connected and displays the appropriate unit of measure. Attach electrode before turning the PH180 on.

## Powering ON/OFF the meter

If the batteries are weak, the '  ' indicator appears on the LCD. Press the ON/OFF key to turn the PH180 on or off. The auto power off feature shuts the PH180 off automatically after 10 minutes of inactivity to preserve battery life.

# Operation

---

## Overview

When the electrode is placed in a solution, the main display and bargraph indicate the pH reading while the lower display reads temperature (readings flash until they have stabilized). The bargraph is 'center zero', i.e. at pH 7 there is no display. As the pH rises, the bar moves from the center to the right. If the pH drops, the bar moves from the center to the left.

## pH Calibration (1, 2, or 3 points)

A two point calibration with a buffer of 7 plus 4 or 10 (whichever is nearest to the expected sample value) is always recommended. A one point calibration (choose the value closest to the expected sample value) is also valid.

1. Place the electrode into a buffer solution (4, 7, or 10). Press and hold the **CAL/RECALL** key until "CAL" appears in the lower (temp.)

display. When doing a 2 or 3 point calibration, calibrate with pH 7 buffer first, then follow with pH 4 then the pH 10 buffer.

2. The WQ120<sup>®</sup> automatically recognizes the solution and calibrates itself to that value (the circled number on the LCD will match the solution). Note that if the solution is more than 1 pH unit off from the L (4), M (7), or H (10) pH buffer, or if the electrode slope is low, the WQ120<sup>®</sup> will assume an error and abort the calibration ('End' will be displayed, and the unit will return to measure mode.)
3. During calibration, the pH reading flashes on the main display.
4. When calibration is complete, the WQ120<sup>®</sup> automatically displays "SA", then "End" and returns to normal operation mode.
5. The appropriate circled indicator (L, M, or H) appears on the LCD when a particular calibration or series of calibrations has been completed within one power on cycle. When the WQ120<sup>®</sup> is turned off, the circled indicator configuration and the calibration data will be retained.
6. For a two or three-point calibration, repeat steps 1-4.
7. See **Reset Calibration Data** to clear all calibration data from the meter.

## Changing the Displayed Temperature Units

To change the displayed temperature units (°C or °F):

1. With the unit OFF, press and hold down the **CAL/RECALL** button.
2. With the **CAL/RECALL** button depressed momentarily press the **ON/OFF** button. When "SELF CAL" appears in the display release the **CAL/RECALL** button. The unit will power on with temperature displayed in the new units.



## Data Hold

Press the **MODE/HOLD** button to hold (freeze) a reading in the display. The meter will enter the HOLD mode and the "HOLD" indicator will appear.

Note: This also stores the reading.

Press the **MODE/HOLD** button again to return to normal operation.

## Storing Readings into Memory


1. Press the **MODE/HOLD** button to store a reading. The storage location number will be displayed on the lower display, while the main display shows the stored reading. The meter will enter the HOLD mode and the "HOLD" indicator will appear.
2. Press the **MODE/HOLD** button again to exit the HOLD mode and return to normal operation.
3. If more than 25 readings are stored, previously stored readings (starting with number 1) will be overwritten.



## Recalling Stored Readings

1. Press the **CAL/RECALL** button and then press the **MODE/HOLD** button. A location number (1 through 25) will briefly appear and then the value stored in that location will appear. The displayed units will flash, indicating that the storage recall mode is active.



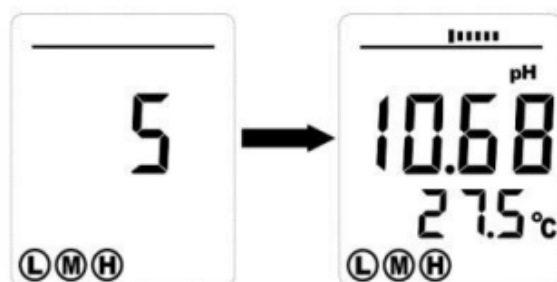

2. The last stored reading will be displayed first. Pressing and releasing the **MODE/HOLD** button will scroll through the stored readings one at a time. The location number is displayed first, followed by the reading stored in that location.
3. To exit the storage mode, press the **CAL/RECALL** button and the meter will return to normal operation, after displaying "End".

## Clear Stored Memory

With the unit on, press and hold ON/OFF for 4 seconds. "CLr" will be briefly displayed when the memory is cleared.

## Auto Power OFF

The auto power off feature automatically shuts the meter off 10 minutes after the most recent button press.

## Auto Power OFF Disable

To disable the Auto Power Off feature:

1. Turn the unit on
2. Press **CAL/RECALL** once (**Quickly**)
3. Immediately and simultaneously press the **MODE/HOLD** and **ON/OFF** buttons for approximately 2 seconds, until "oFF" is briefly displayed

To disengage this feature, turn the unit off with the **ON/OFF** button. The next time the unit is powered up, Auto Power OFF mode will be engaged again.

## CAL Reminder Display

When the PH180 is turned on in the pH mode for the 15th time without recalibration, the 'CAL' icon appears on the LCD indicating that the PH180 may require calibration. Some applications may require recalibration of the electrode more frequently than others. The CAL display is simply a reminder and will turn off when the pH electrode is recalibrated.

## RENEW Display

A flashing 'RENEW' warning indicates that the probe is not performing to expected specifications. If cleaning and recalibration does not cause the RENEW icon to disappear, replace the probe (see optional accessories on the last page of this manual). The RENEW display appears as a result of the pH electrode slope falling below 65% of a nominal slope.

## Measurement and Display Considerations

If the unit appears to be locked (display frozen). It is possible that the Data Hold mode has been inadvertently accessed by pressing the **MODE/HOLD** button. ("HOLD" will be displayed in the bottom left of the LCD.) Simply press the **MODE/HOLD** button again or turn the meter off and then on.

For maximum accuracy, allow sufficient time for the temperature of the probe to reach the temperature of the sample before calibrating. This will be indicated by a stable temperature reading on the display.

## Reset Calibration Data

Follow this procedure to clear all calibration data from the meter. Resetting the calibration data may be necessary when new calibration solutions are used or accuracy of measurements is in question.

1. Turn off the meter.
2. Press and Hold the Cal/Recall and Mode/Hold buttons.
3. Momentarily press the On/Off button, as soon as the display comes on, release all 3 buttons.
4. The display will show "dFLt rSt" (default reset) and all of the calibration data will be erased. If "dFLt rSt" does not appear, retry the procedure.

Proceed to the calibration routine for pH and Conductivity.

## Battery Replacement

---

1. Twist off the battery compartment cap
2. Replace the four (4) 2032 batteries observing polarity.
3. Replace the battery compartment cap



Never dispose of used batteries or rechargeable batteries in household waste.

As consumers, users are legally required to take used batteries to appropriate collection sites, the retail store where the batteries were purchased, or wherever batteries are sold.

**Disposal:** Do not dispose of this instrument in household waste. The user is obligated to take end-of-life devices to a designated collection point for the disposal of electrical and electronic equipment.


## Other Battery Safety Reminders

- Never dispose of batteries in a fire. Batteries may explode or leak.
- Never mix battery types. Always install new batteries of the same type.



## Specifications

---

Display	Multifunction LCD with Bargraph
Operating conditions	32 to 122°F (0 to 50°C) / < 80% RH
Range and Accuracy	0.00 to 14.00 / $\pm 0.01$ pH typical
Temp. Compensation	Automatic from 32 to 194°F (0 to 90°C)
Temperature Range	32 to 194°F (0 to 90°C)
Temperature Resolution	0.1° up to 99.9 then 1° thereafter
Temperature Accuracy	$\pm 1^{\circ}\text{C}/1.8^{\circ}\text{F}$
Measurement storage	25 tagged (numbered) readings
Power	Four (4) CR2032 button batteries
Low battery indication	'  ' appears on the LCD
Auto power off	After 10 minutes of inactivity
Dimensions	1.5 x 7.8 x 1.5" (38 x 198 x 38 mm)
Weight	3.5 oz (100 g)

## Warranty

---

Triplett / Jewell Instruments extends the following warranty to the original purchaser of these goods for use. Triplett warrants to the original purchaser for use that the products sold by it will be free from defects in workmanship and material for a period of (1) one year from the date of purchase. This warranty does not apply to any of our products which have been repaired or altered by unauthorized persons in any way or purchased from unauthorized distributors so as, in our sole judgment, to injure their stability or reliability, or which have been subject to misuse, abuse, misapplication, negligence, accident or which have had the serial numbers altered, defaced, or removed. Accessories, including batteries are not covered by this warranty

**Copyright © 2021 Triplett**  
**[www.triplett.com](http://www.triplett.com)**

***DIREKTRONIK***

**Kontakta Sverige:**  
[www.direktronik.se](http://www.direktronik.se)  
[info@direktronik.se](mailto:info@direktronik.se)  
Ring: 08 52 400 700

**Kontakt Norge:**  
[www.direktronik.no](http://www.direktronik.no)  
[info@direktronik.no](mailto:info@direktronik.no)  
Ring: 55 50 91 50